



X-pendable Cartridge USE, DURATION, & STORAGE

GENERAL

Like a magnet, the purpose of our X-pendable cartridges is to attract and capture various contaminants.

This is facilitated through an assortment of sorbents. The ultimate goal is to capture as much of these contaminants as possible.

But beware, their capacity is limited. Leave them exposed to the atmosphere and they will do their job ...until full and their lifespan can easily shrink prematurely. Your goal with these cartridges is to **protect them until ready for use.**



DURATION

Useable duration (aka, Processing Capacity) is affected by variables such as:

- ❖ The type of process gas
- ❖ Temperature, pressure, and nature/level of contaminants
- ❖ Types and volume of sorbents
- ❖ Dwell time

Minor changes in any of these could drastically impact the overall capacity and useable lifespan of the cartridge(s). Digital monitor can aid in watching the duration of the cartridge.

(LabOnLocale is such a device

www.LabOnLocale.com)

Maximum usable life is 6 months once installed

STORAGE

X-pendable's carry a guaranteed 2-year shelf life in our standard package, so long as unopened.

We also offer a 4-year package option. The manufacture date is encoded in the Lot Number. Contact us for the exact manufacture date.

USE

See instructions included with your cartridge. Change all cartridges in system at same time. Not for use with Oxygen unless noted. Change cartridges once LifeBand has turned beige/pink



Others make air-&-gas...

We ensure that it's PURE

4790 NW 157th Street, Miami FL 33014 USA
305-430-0550 info@Lawrence-factor.com
www.Lawrence-Factor.com